


防腐剤の分析-2

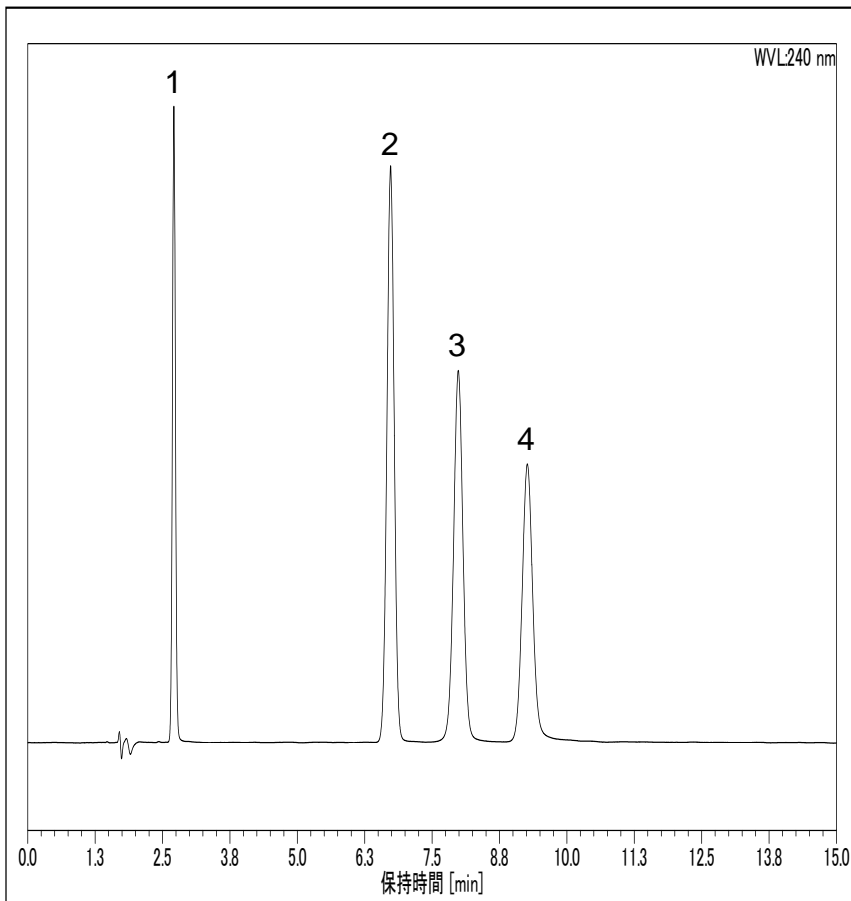

Sample

1. *p*-Hydroxybenzoic acid
(0.11mg/ml)

2. Benzoic acid (0.30mg/ml)

3. Salicylic acid (0.24mg/ml)

4. Dehydroacetic acid (0.61mg/ml)

Conditions;

Column: Develosil XG-C18M (5 μ m) 4.6x150mm

Mobile phase: Acetonitrile/0.2%Formic acid=25/75

Flow rate: 1.0ml/min

Temperature: 40°C

Detection: UV240nm

Injection volume: 2.0 μ L

System: DIONEX/Ultimate 3000 Series